



Armin Baack

Bedarf und Technik für Labor und Medizin

Ihr kompetenter Partner für Ihr Labor

Thermotrack Mobile Edition

MAIN FUNCTIONS

- Read Thermo and Hygro Buttons on your Android Smartphone or Tablet
- Verify alarms
- Relaunch Mission

Advantages

- On site mobile controls
- Easy to use
- Synchronize with other Thermotrack applications or your server
- Free download

Description

Thermotrack Mobile Edition is a free app which enable you to use your Thermo and Hygro Buttons anywhere using your Android Smartphone or Tablet.

Thermotrack Mobile can:

- Control occurrence of Temperature alarms on site
- Download data on your smartphone
- Transfer it on your favorite application
- Relaunch the current mission of the Thermo Button or Hygro Button.

Download the application from the Google Play*. It's free! You just need a Thermo Button or Hygro Button reader which you can get from your PlugAndTrack reseller.



Alarm CONTROL

Plug the reader on a Thermo Button or Hygro Button and touch the icon «alarm». The icon turns:

Green if no alarms

Orange if temperatures threshold have been reached but within an acceptable duration (Temporization)

Red if both thresholds and temporization have been reached

Workswith

Thermotrack PC
Thermotrack-Online
Thermotrack-Webserve



* Check compliance of your device from the Google Play Store

Download and transfer

Data can be downloaded on your smartphone and transferred to your favorite temperature tracking application, so you can analyze, print and keep it. You also can get them as XML and integrate it into any other software.

If you don't have any GSM connection, then use the WiFi when you're back in your office

mission RELAUNCH

If you want to clear the Button memory and start a new mission, just press the icon «Relaunch»